

Report Generator Options available (Report Properties):

The screenshot shows the 'Report Options' dialog box with the 'General' tab selected. The dialog has a blue title bar with a close button. Below the title bar are five tabs: 'General', 'Filters', 'Stability Summary', 'Stability Summary (cont.)', and 'Alternate Stability Summary'. The 'Options' section contains two text input fields: 'Company Name' and 'Alternate Title'. At the bottom are 'OK' and 'Cancel' buttons.

Report Options

General | Filters | Stability Summary | Stability Summary (cont.) | Alternate Stability Summary

Options

Company Name

Alternate Title

OK Cancel

The screenshot shows the 'Report Options' dialog box with the 'Filters' tab selected. The dialog has a blue title bar with a close button. Below the title bar are five tabs: 'General', 'Filters', 'Stability Summary', 'Stability Summary (cont.)', and 'Alternate Stability Summary'. The 'Storage Condition' section has three checked items: 'not stored', '25°C ± 2°C/60% RH ± 5% RH', and '40°C ± 2°C/75% RH ± 5% RH'. The 'Test Categories' section has four checked items: 'Assay and Related Substances', 'Physical Examination', 'Dissolution', and 'Hardness'. The 'Intervals' section has six checked items: 'Initial', '2 week', '1 month', '3 month', '6 month', and '12 month'. There are scroll arrows on the right side of the 'Intervals' list. At the bottom are 'OK' and 'Cancel' buttons.

Report Options

General | Filters | Stability Summary | Stability Summary (cont.) | Alternate Stability Summary

Storage Condition

- not stored
- 25°C ± 2°C/60% RH ± 5% RH
- 40°C ± 2°C/75% RH ± 5% RH

Test Categories

- Assay and Related Substances
- Physical Examination
- Dissolution
- Hardness

Intervals

- Initial
- 2 week
- 1 month
- 3 month
- 6 month
- 12 month

OK Cancel

Report Options [X]

General | Filters | **Stability Summary** | Stability Summary (cont.) | Alternate Stability Summary

Options

Show Approved Only
 Show Units After Results
 Add Signature Section
 Show Shading
 Show Comments
 Add Notebook Reference
 Show Abbreviated Header
 Show LWR information
 Show Scheduled Status
 Restrict Comments
 Special Instructions
 Show Alternate Title
 Group Packaging Items
 Add Company Name

Show Pull Dates:

Show Test Dates:

Number of Intervals Per Page:

OK Cancel

Report Options [X]

General | Filters | Stability Summary | **Stability Summary (cont.)** | Alternate Stability Summary

Header Options

Left Column	Right Column
<input checked="" type="checkbox"/> Show Product/Strength	<input checked="" type="checkbox"/> Show Lot #/ Mfg. Date
<input checked="" type="checkbox"/> Show Mfg. By/Site	<input checked="" type="checkbox"/> Show Packager/Site
<input checked="" type="checkbox"/> Show Batch Size	<input checked="" type="checkbox"/> Show Packaging Date
<input checked="" type="checkbox"/> Show Purpose of Study	<input checked="" type="checkbox"/> Show Study Start Date
<input checked="" type="checkbox"/> Show Expiration Date	<input checked="" type="checkbox"/> Show Study Duration
<input checked="" type="checkbox"/> Show Drug Substance Information	<input checked="" type="checkbox"/> Show Package Item Information
<input type="checkbox"/> Show Test Locations	

OK Cancel